

## **DAPI Staining**

### **Stock Solution:**

5  $\mu$ L DAPI (10 mg/mL solution)

995  $\mu$ L PBS

\*\*store at 4°C in the dark

\*\*solution is viable for 4 months

### **Working solution:**

\*\*prepare when ready to use

10  $\mu$ L Solution 1

990  $\mu$ L PBS

- Bake slides overnight at 40°C
- Incubate slides with working solution for 1-5 minutes
- Wash 2 x for 10 minutes each with PBS
- Coverslip slides with Mowoil